



CAL4-E

User Manual

5. May 2017





Content

1 INTRODUCTION	3
1.1 Conformity	3
1.2 Intended use.....	3
1.3 General safety information	3
2 GETTING STARTED	3
2.1 Scope of delivery	3
3 OPERATING INSTRUCTIONS HARDWARE	3
3.1 Requirements	3
3.2 Connection to the projector	4
3.3 Starting the system.....	4
3.4 Endoscopy positioning	5
4 ADDITIONAL INFORMATION	5
4.1 Care instructions	5
4.2 Disposal instructions	5
5 TECHNICAL DATA SHEET	5



1 INTRODUCTION

Important information: Read the manual carefully before using this device.

Inappropriate utilization may cause damages to the device, to the DUT (device under test) and/or other components of your setup.

Keep these instructions in a safe place and pass them to any future user.

1.1 Conformity

We, Image Engineering GmbH & Co. KG, hereby declare, that the CAL4-E corresponds to the essential requirements of the following EC directive in its current version:

- Electromagnetic Compatibility - 2014/30/EU

1.2 Intended use

The integrating sphere is designed for the measurement of color, resolution, OECF, dynamic range and noise when using the endoscopy light source.

- Only suitable for indoor use.
- Place your system in a dry and constant tempered environment without any interfering light.
- The optimal ambient temperature range is 22 to 26 degrees Celsius.

1.3 General safety information

- **Do not** look directly in the open sphere or light source when using high intensities.
- **Do not** open the device without any prior instructions from the Image Engineering support.

2 GETTING STARTED

2.1 Scope of delivery

CAL4-E - 30cm integrating sphere (without light source)

Four adapters for different kind of endoscopes

High temperature resistant cold-light cable, XENON approved

3 OPERATING INSTRUCTIONS HARDWARE

3.1 Requirements

- Endoscope
- Projector



3.2 Connection to the projector

Connect the CAL4-E to your projector by using one of the four adapters with the fiber.



3.3 Starting the system

Use the brackets to place a test charts in front of the CAL4-E orifice and turn on the projector's light source.



NOTICE

The CAL4-E device can only operate with high precision, when the projector's light source is ready to use. Please consult the user manual of your projector to get the information about the warmup time for a constant illuminance.



3.4 Endoscopy positioning

Please verify that

- the image height includes the test chart height.
- the lens is exactly in the middle of the test chart

Not fulfilling these requirements will lead to an inhomogeneous illuminated field of view, or might provide questionable measurement results.

4 ADDITIONAL INFORMATION

4.1 Care instructions

- Do not touch, scratch or pollute the diffuser.
- If there is any dust on the diffuser clean it with an air blower.
- Do not remove the fiber from the CAL4-E, otherwise the illumination is invalid
- Store and transport the CAL4-E only in the delivered hard case.

4.2 Disposal instructions

After the service life of CAL4-E, it must be disposed properly. Observe your national regulations. Make sure that CAL4-E cannot be used by third parties after disposing of it.

Contact Image Engineering if assistance for disposal is required.

5 TECHNICAL DATA SHEET

See annex for the technical data sheet. It can also be downloaded from the website of Image Engineering : www.image-engineering.com.